


New Standards in Fibrosis Quantitation: The Value of Label-Free Second Harmonic Generation Imaging

Li Chen and Mathieu Petitjean

Genesis Imaging Services, PharmaNest, Princeton, NJ, USA

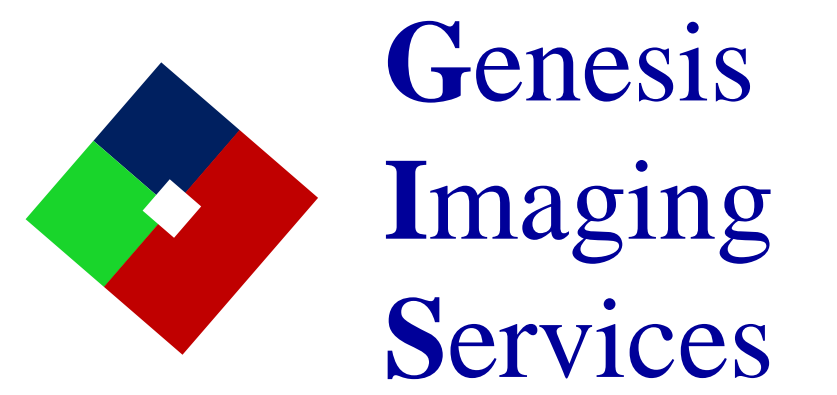

Background


Fibrosis is associated with excessive accumulation of extracellular matrix in response to persistent injury, inflammation, and abnormal wound healing. The current standard to assess fibrosis are conventional staining and histopathological criteria scoring. This assessment has limitation as it uses a narrow range scoring (when scoring exists), it is prone to observer variations, and requires multiple histopathology workflows and stains.

- Here, we use second harmonic generation (SHG) together with two-photon excitation fluorescence (2PE) imaging and computerized image analysis algorithms to provide a novel, sensitive, and efficient method for collagen quantitation. We extract information on the collagen fibers in various animal models to assess fibrosis as it advances or regresses in response to therapeutic compounds.
- We present animal models including NASH livers, IPF lungs and UVO kidneys treated with or without reference anti-fibrotic drugs including Nintedanib and Pirfenidone.
- In addition to basic quantifiable metrics including total collagen, we offer novel analysis capabilities that conventional method are not able to provide such as tissue regional segmentation (for livers, lung, kidney), collagen network structure, collagen fiber density, etc. These metrics maybe the key to better fibrosis scoring and staging.
- Due to the nature of the technology, this stain-free SHG/2PE imaging and automated image analysis allow for a reliable and quick turnaround time for data results.

*We thank GenScript and Intercept Pharmaceuticals for providing the conventional staining images and respective graphs. We thank HistoIndex and A*STAR (Singapore) for their assistance in the image analysis of the IPF rodent model images.

METHOD

Tissue Preparation, Instrumentation, and Workflow

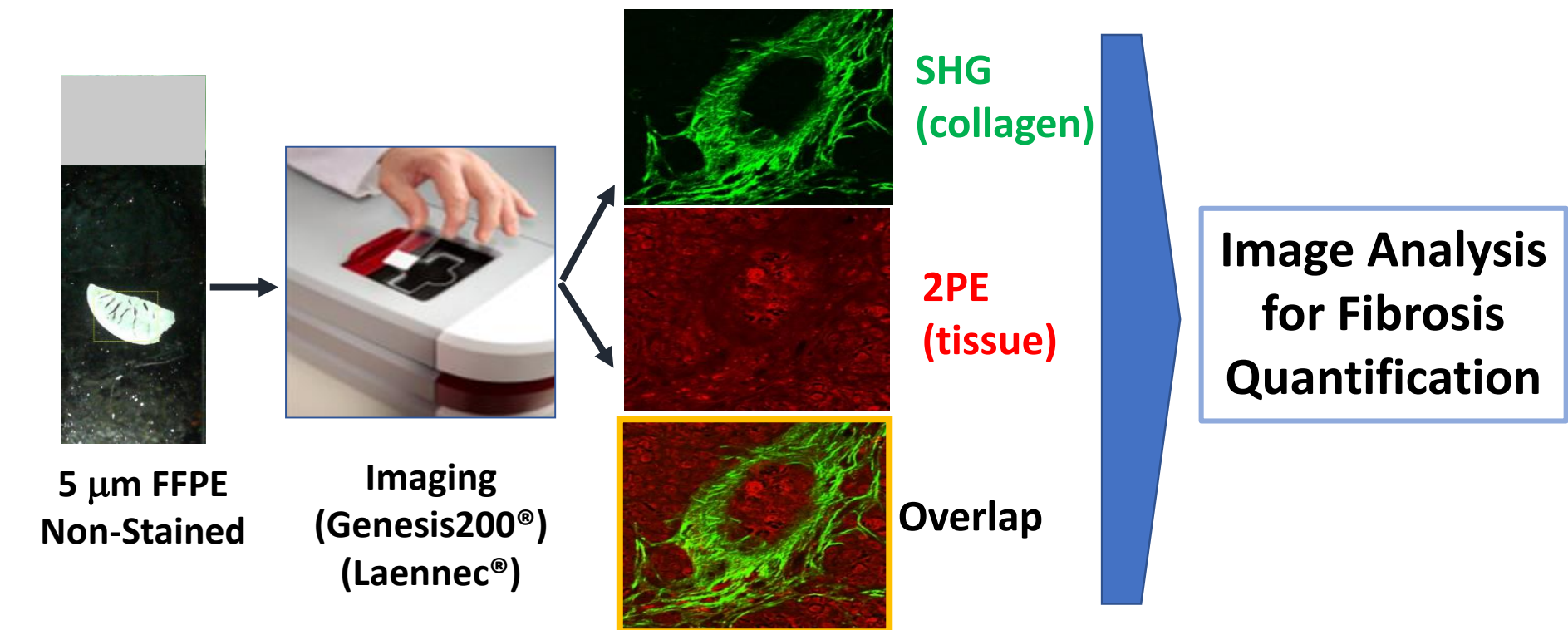

- 5-200μm FFPE or Frozen sections
- Stain-Free and label-free imaging
- Fully quantitation of collagen fibers
- High Resolution (0.39um @ 20X)
- Non-destructive (tissue re-usable)
- 2-Photon Excitation (2PE): auto fluorescence to delineate tissue morphology and injury
- Second Harmonic Generation (SHG): Collagens I and III
- Image Analysis Software optimized and validated with pathologists

Amylin Liver NASH Model

Leptin-deficient (*lep^{ob/ob}*) mice were allowed ad libitum access to normal chow (low-fat diet w no fructose nor cholesterol) or to modified ALIOS diet (high trans fat (40%), fructose (22%), cholesterol (2%) in food pellets) with or without therapeutic drugs.

IPF Lung Model

C57BL/6 mice were subjected to intratracheal instillation of bleomycin for induction of lung epithelial injury, inflammation, and fibrosis. Pirfenidone or Nintedanib was given 7days after start of bleomycin (21days study).

UVO Kidney Model



SD rats were subject to surgical ligation of ureter of left kidney, while leaving the right contralateral kidney as control. This induces progressive epithelial injury, tubulointerstitial inflammation and fibrosis of the left kidney. Pirfenidone is given 7 days after start of ureter unilateral obstruction (UVO) (28 days study).

Models	Group 1	Group 2	Group 3	Group 4
NASH Livers	Chow	NASH	NASH + OCA	NASH + Drug A
IPF Lungs	Saline	Bleomycin (BLM)	BLM + Nintedanib (NTN)	BLM+ Pirfenidone (PFD)
UVO Kidney	Sham	UVO	UVO+PFD	

- Collagen Area Ratio (CAR): % collagen area within area of interest
- Collagen Reticulation Index (CRI): measures complexity of the collagen fiber network
- Collagen Fiber Density (CFD): collagen density within fiber based on pixel intensity
- Fat area Ratio (FAR): % fat area within area of interest



Figure 1. Modified ALIOS diet-feed mice induces hepatic fibrosis and steatosis treated with or without obeticholic acid (OCA) or Drug B.

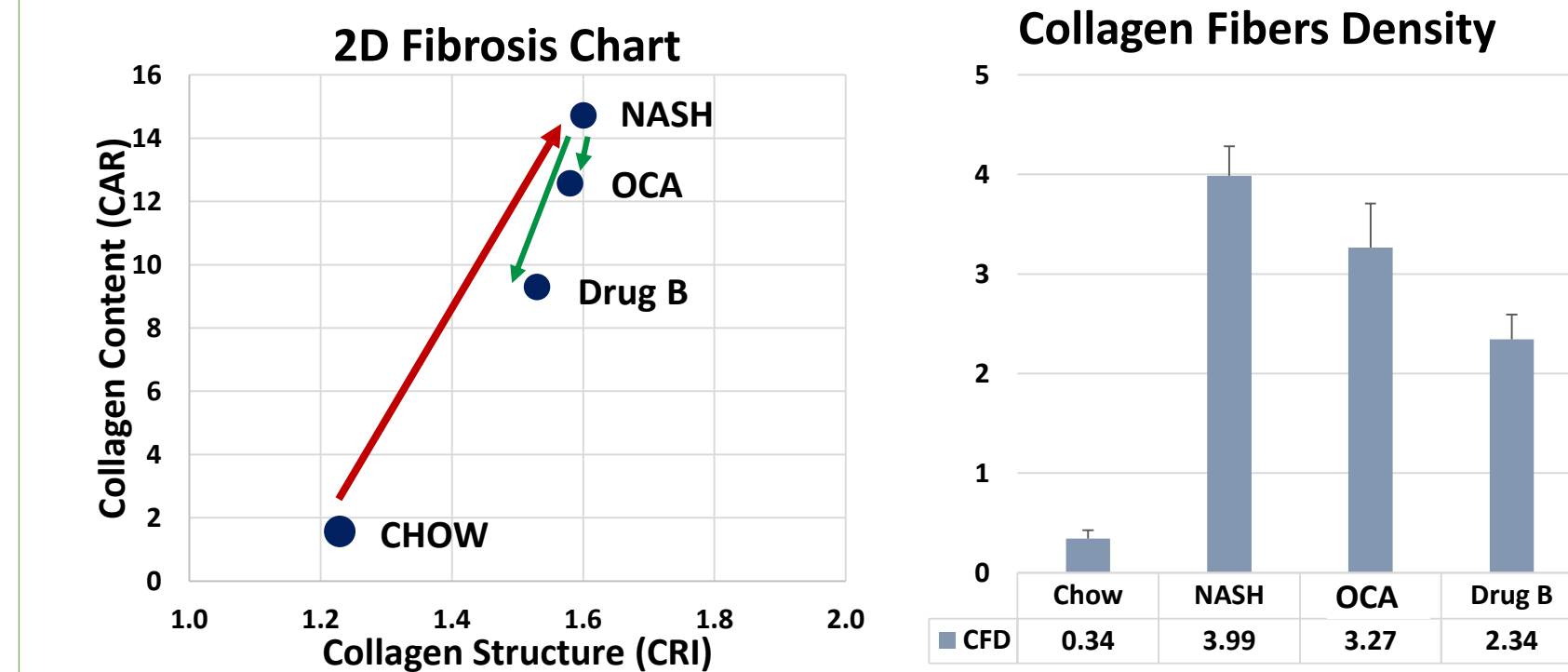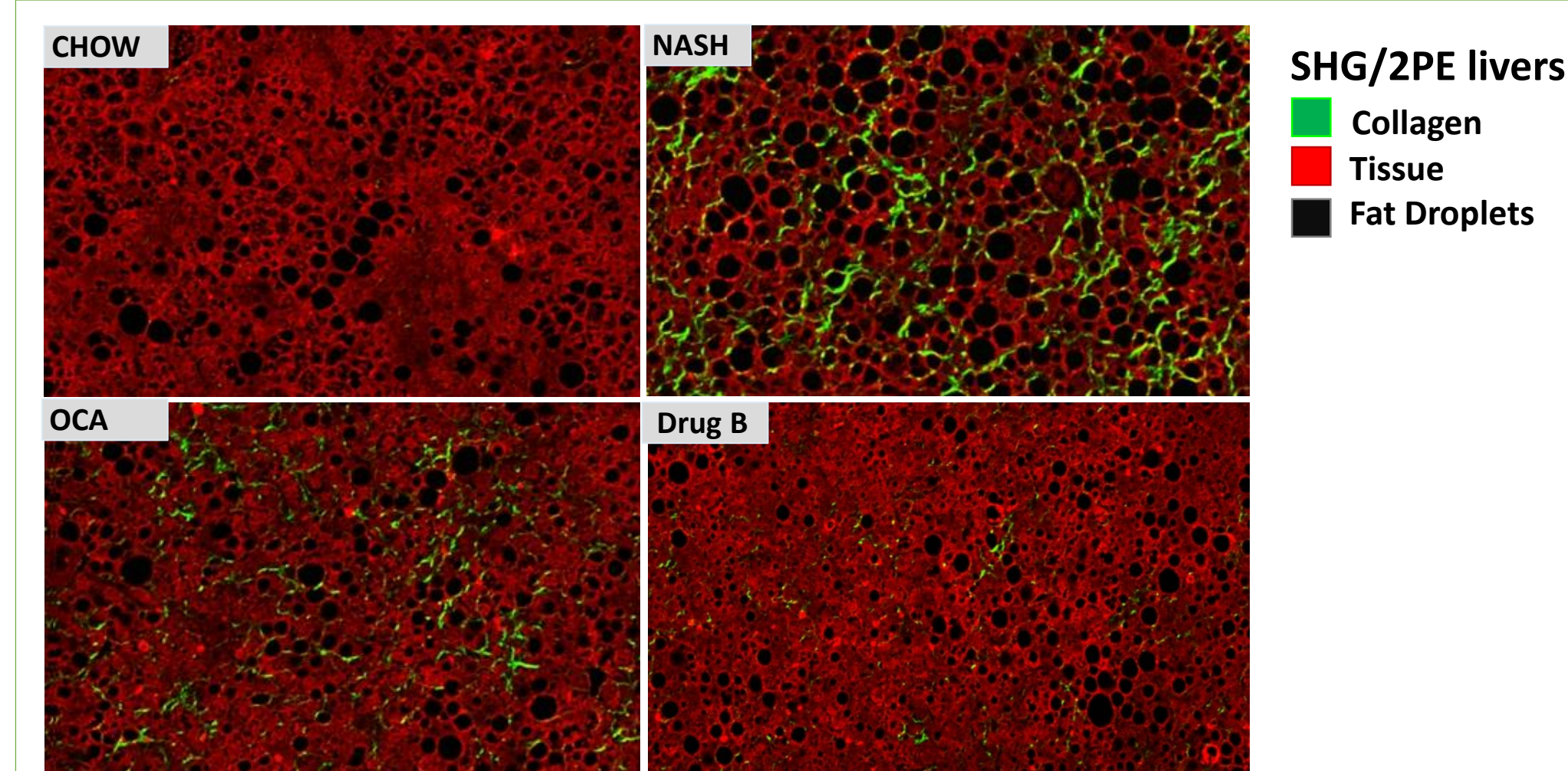

Automated NASH liver image analysis for concurrent fibrosis and steatosis quantitation.



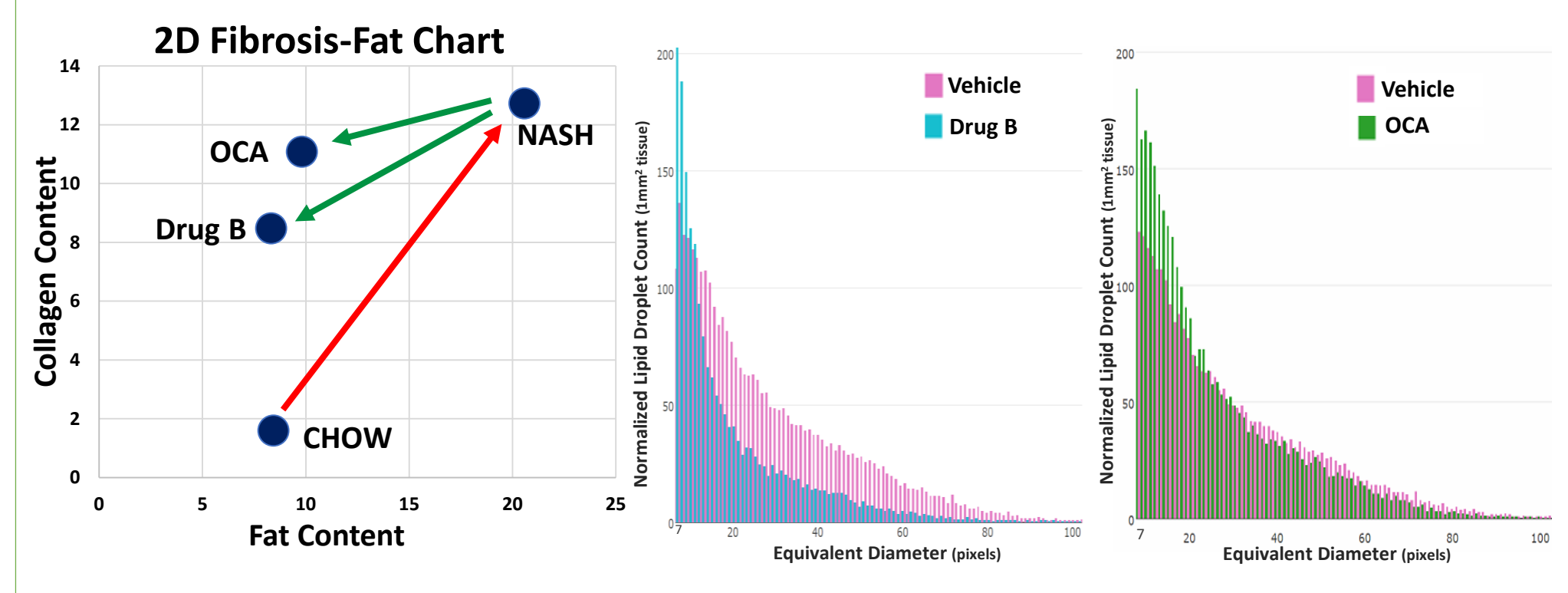


Figure 2. Modified ALIOS diet-feed mice induces steatosis and fibrosis. Conventional staining for comparison. Collagen la1(brown)

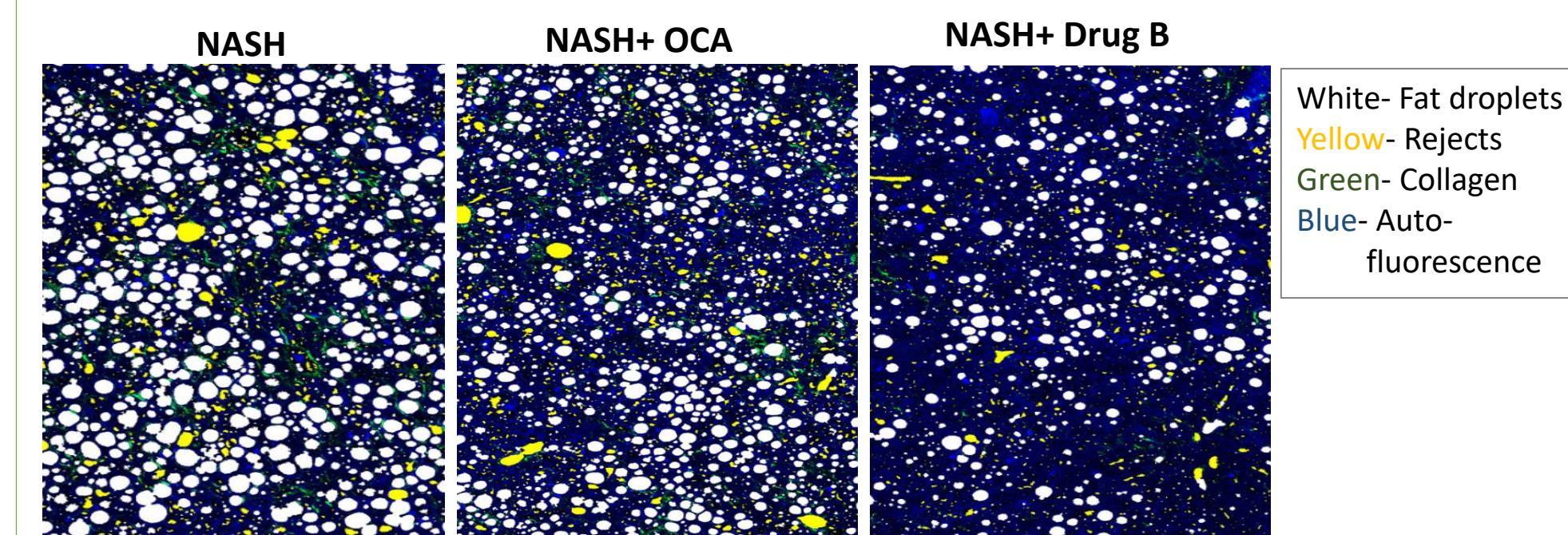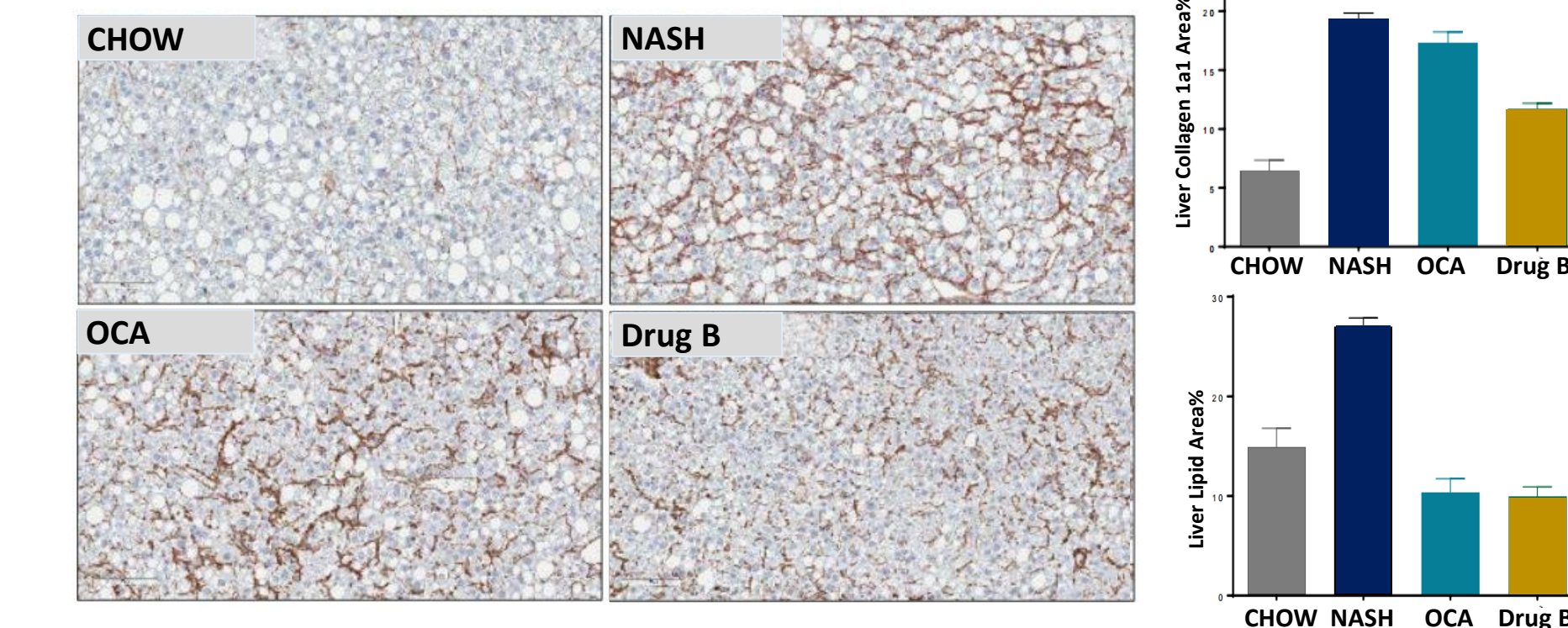

RESULT

Figure 4. Bleomycin induced lung fibrosis and treated with or without Nintedanib or Pirfenidone

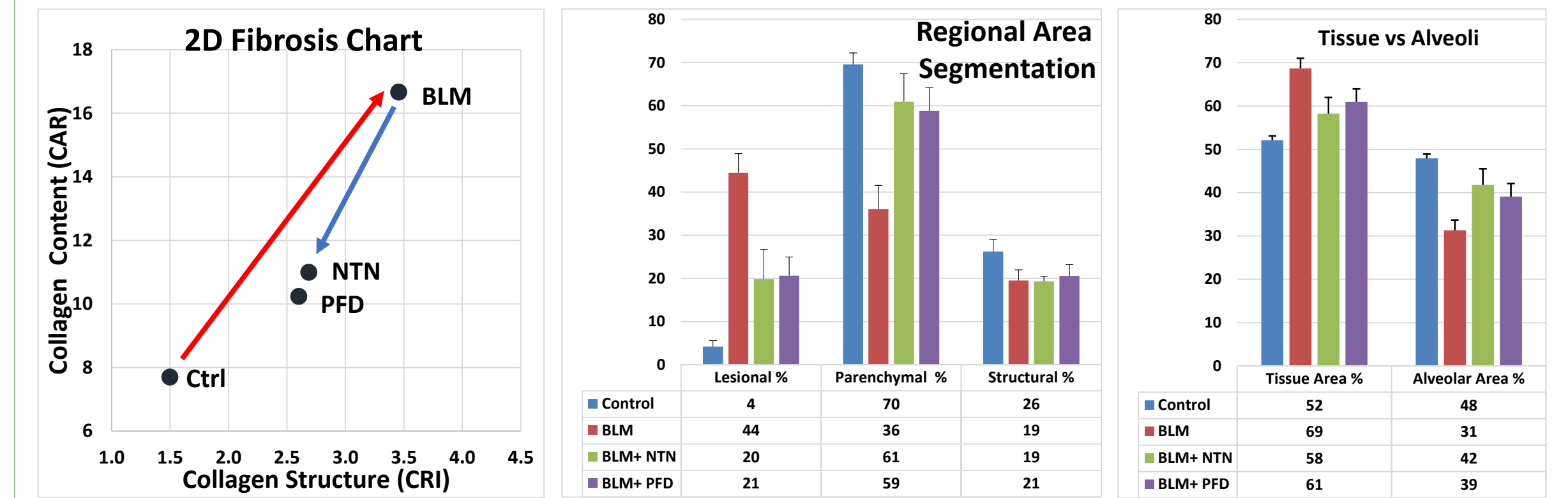
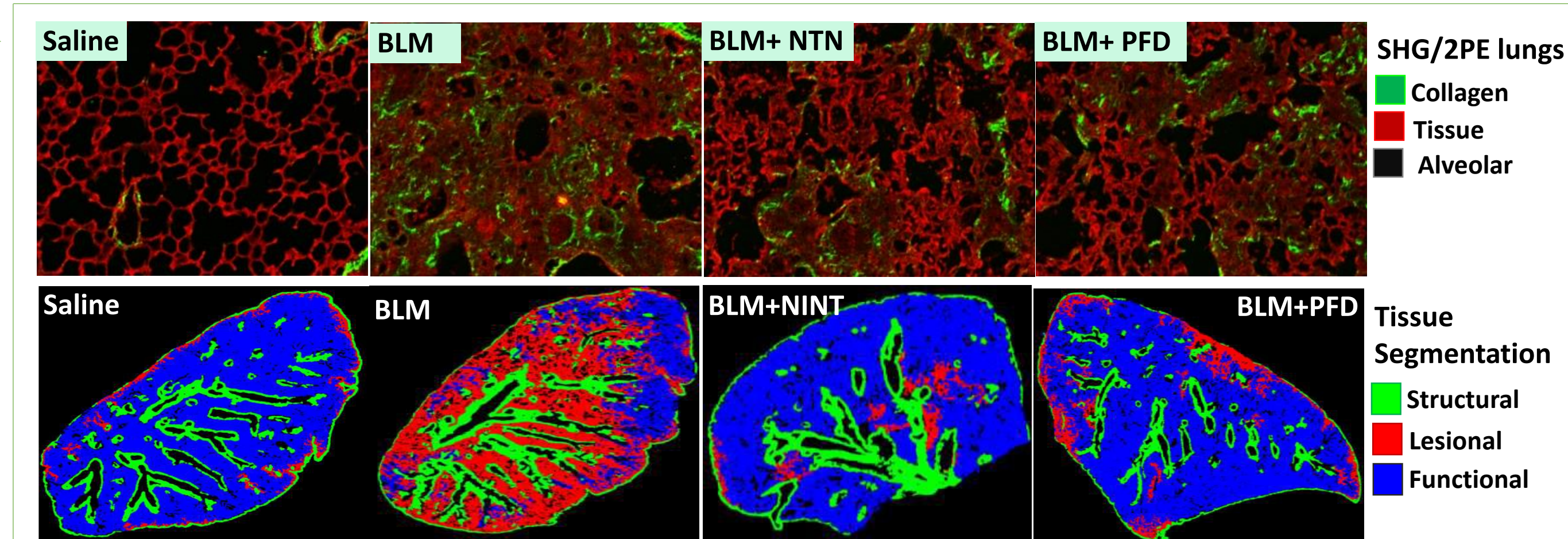


Figure 4. Bleomycin induced lung fibrosis and treated with or without Nintedanib or Pirfenidone.

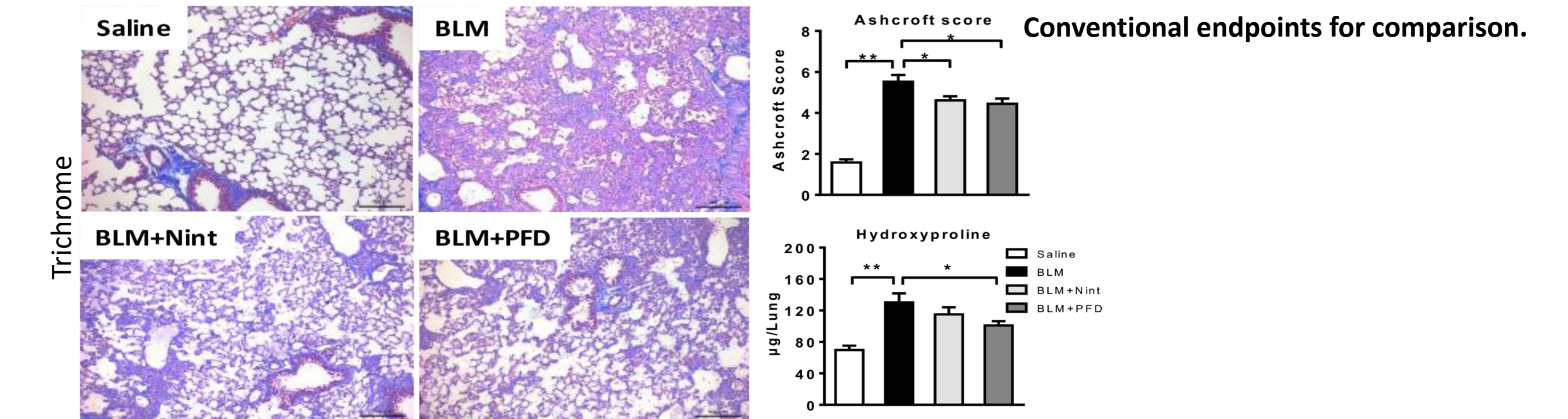


Figure 5. UVO induced kidney fibrosis in rat treated with or without Pirfenidone

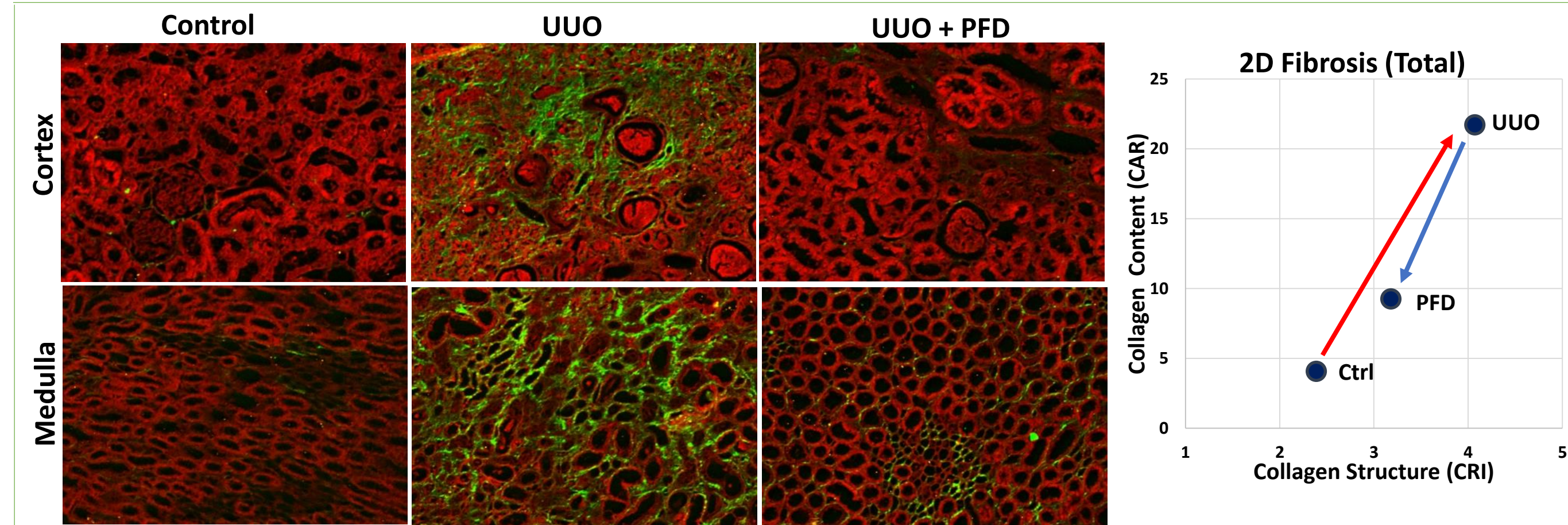


Figure 6. UVO induced kidney fibrosis in rat treated with or without Pirfenidone.

